
7694 Oven Top

	Description	HP Part No.	Qty
1	Spring lock	301-077-HSP	1
2	Spacer	300-520-HSP	2
3	Screw, cover lock	301-147-HSP	2
4	Frame, cover oven	301-112-HSP	1
5	Hinge, oven case	301-119-HSP	1
6	Rod, lever lock cover	301-019-HSP	1
7	Oven top cover	301-111-HSP	1
8	Carousel assembly (See figure 6-8)		1
9	Frame, base case	301-117-HSP	1
Associated screws, nuts, and washers			
A	Screw, M3 X 10 pan head	0515-0925	2
B	Washer, toothed D3	2190-0638	4
C	Washer, flat D3	3050-0891	4
D	Nut, M30535-0004	2	
E	Washer, flat D4	3050-0893	3
F	Washer, toothed D4	2190-0640	3
G	Nut, M40535-0006	3	
H	Screw, socket head M4 X 12	300-837-HSP	3
I	Washer, toothed M4	2190-0640	3
J	Washer, flat M4	3050-0893	3

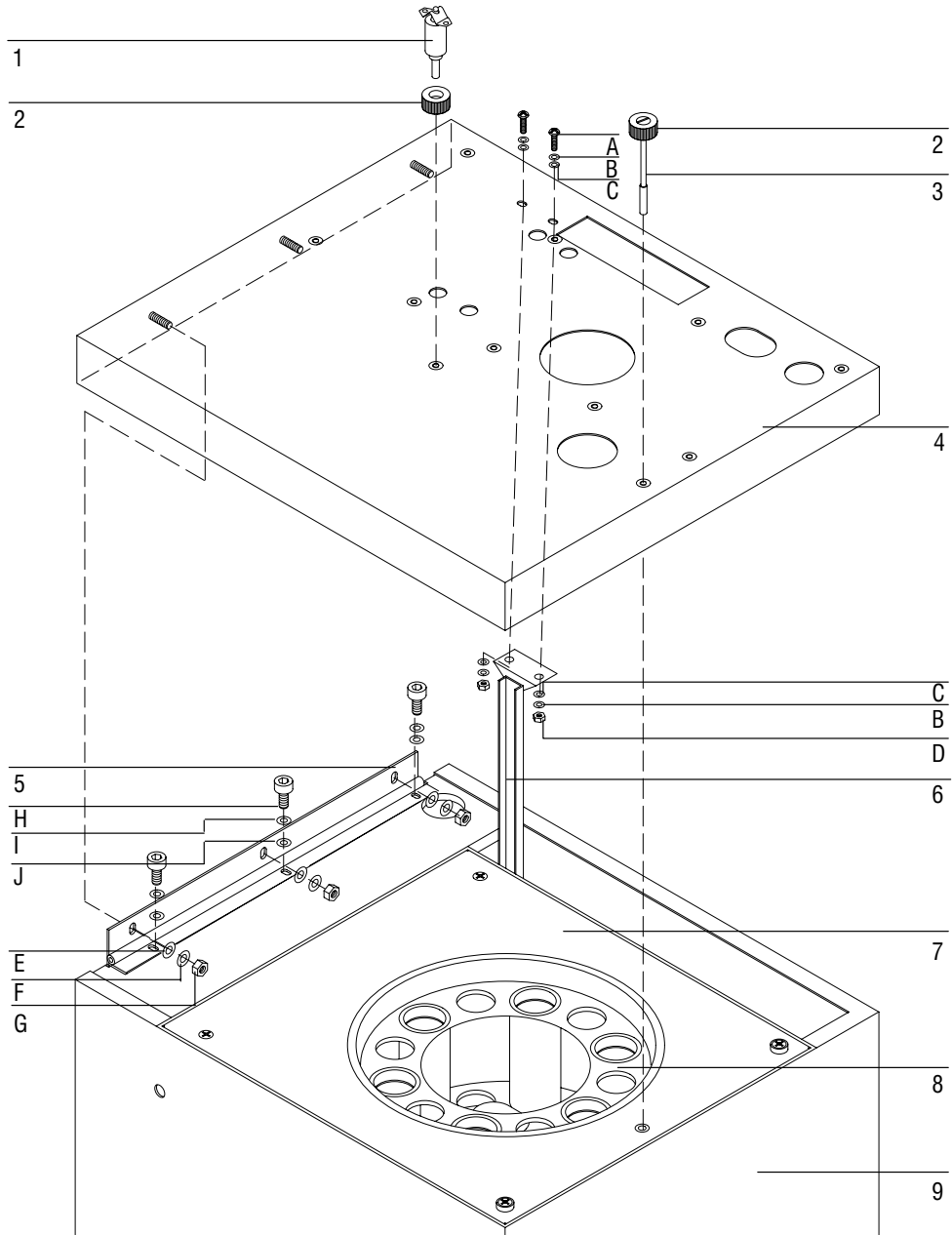


Figure 6-1. Removing the Oven Top