

Chromatography Solutions

Avantor[®] ACE[®] Excel[®] Glycan

UHPLC/HPLC columns for reproducible
glycan analysis



Avantor[®] ACE[®]

GLYCAN ANALYSIS

- Analysis of the glycosylation pattern of biopharmaceutical compounds is essential for their complete characterisation
- Monoclonal antibodies (mAbs) are one of the largest growing areas of biotherapeutic interest and are heavily glycosylated molecules with each mAb containing at least two glycosylation sites as shown in Figure 1 below
- Changes to the glycosylation of a biopharmaceutical can affect the efficacy and safety of the drug, so it is extremely important that it is thoroughly monitored
- Due to the highly polar nature of released glycans, a hydrophilic stationary phase is ideal for their separation
- Labelling with a fluorophore such as 2-aminobenzamide (2-AB) allows the released glycans to be analysed using fluorescence detection

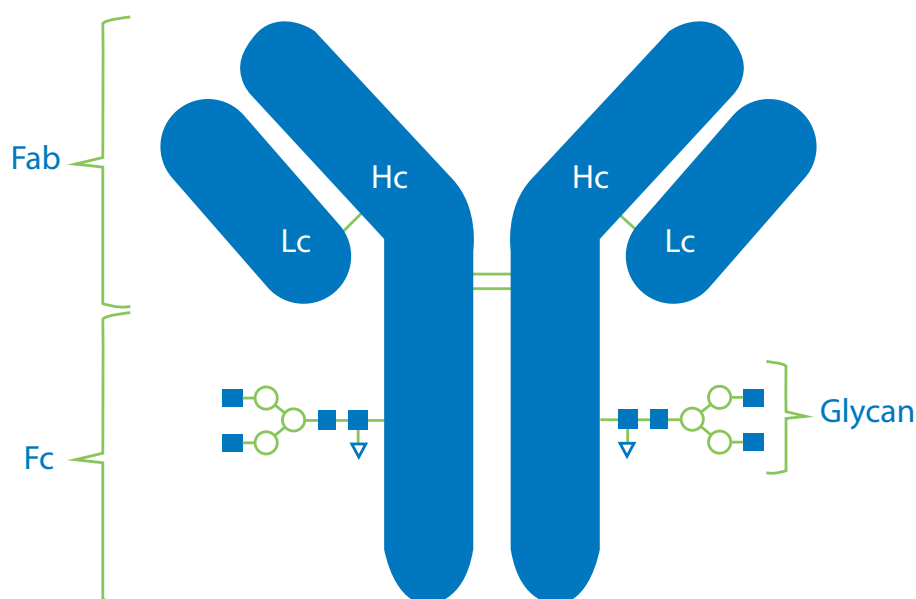


Figure 1: Schematic showing glycosylation regions of a monoclonal antibody.

Phase	Endcapped	Particle size (µm)	Pore size (Å)	Surface area (m ² /g)	Carbon load (%)	Recommended pH range
ACE® Excel® Glycan	No	1.7, 3	100	300	7.0	2.0 - 7.0

To extend column lifetime ACE UHPLC guard cartridges (Cat. No. EXL-1P16-GD2U / Cat. No. EXL-1P16-GD4U) are recommended.

For the guard system, guard cartridge holder (H0011) is also required.


For HPLC column connections up to 6000 psi, PEEK™ fingertight fittings (Cat. No. ACE-CC10) are recommended - for further details please contact chromsupport@avantorsciences.com

For UHPLC column connections up to 25000 psi, reuseable fittings (Cat. No. EXL-CC10) are recommended - for further details please contact chromsupport@avantorsciences.com or visit vwr.com/ace

GUARANTEED REPRODUCIBILITY

- The ACE® Excel® Glycan column provides excellent batch-to-batch reproducibility as shown below
- The dextran ladder standard consists of a mixture of glucose homopolymers of between 1 to 15 monomer units, and is a useful tool in identifying more complex glycoforms based upon correlated retention times

BATCH-TO-BATCH REPRODUCIBILITY


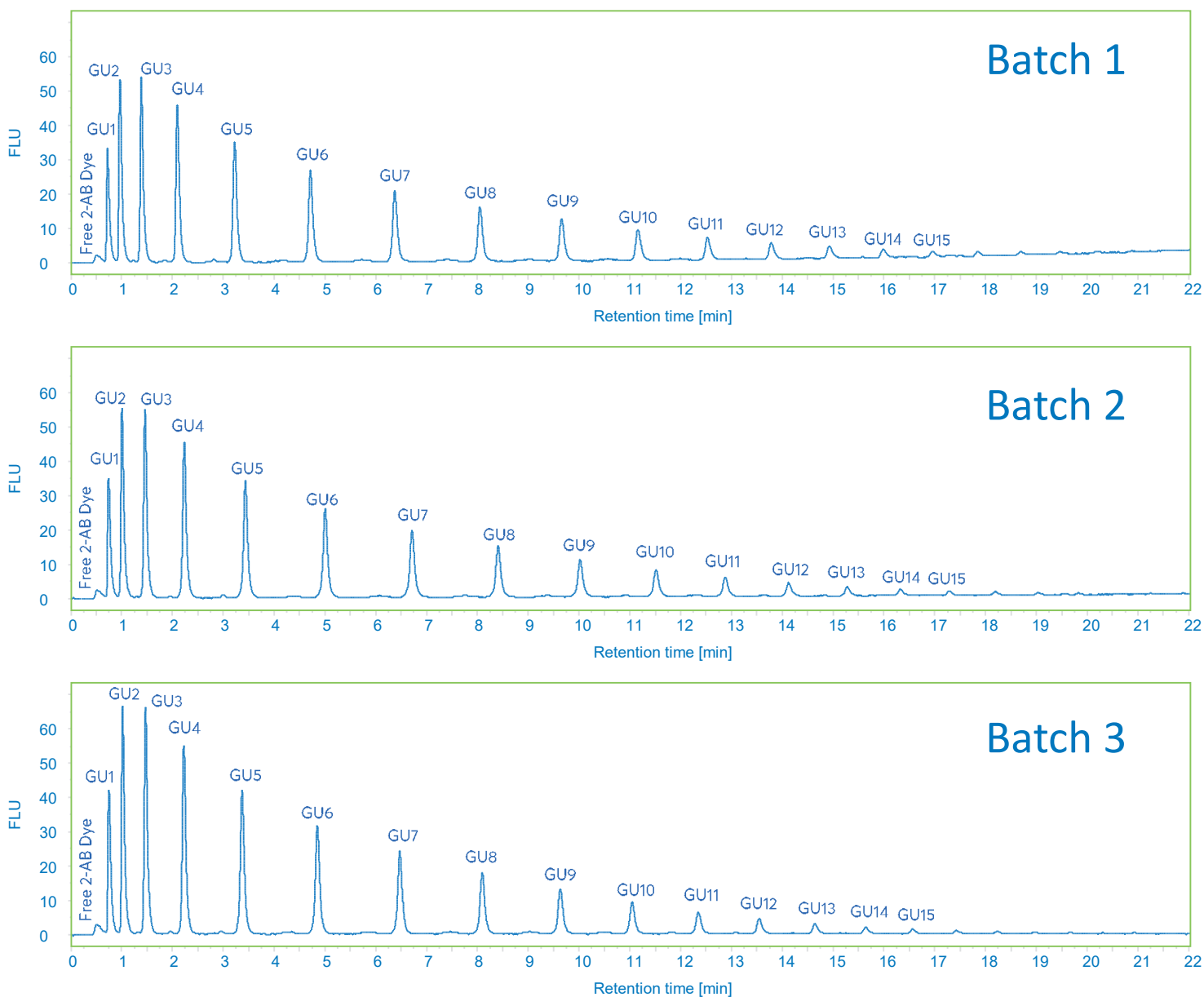


Sample: 2-AB labelled Dextran Ladder standard
Column: ACE Excel 1.7 Glycan, 100x2.1 mm
Cat. No.: EXL-1716-1002
Mobile phase: A = 100 mM ammonium formate (pH 4.5) and B = MeCN
Gradient: 75 – 52.5% B in 24 min
Flow rate: 0.5 ml/min
Temperature: 55 °C
Injection: 1 µl
Detection: FLD, λ_{ex} = 260 nm, λ_{em} = 430 nm

GUARANTEED SCALABILITY

The ACE® Excel® Glycan column provides exceptionally reproducible scalability between the 1.7 μm and 3 μm particle sizes, as demonstrated below

PARTICLE SIZE COMPARISON

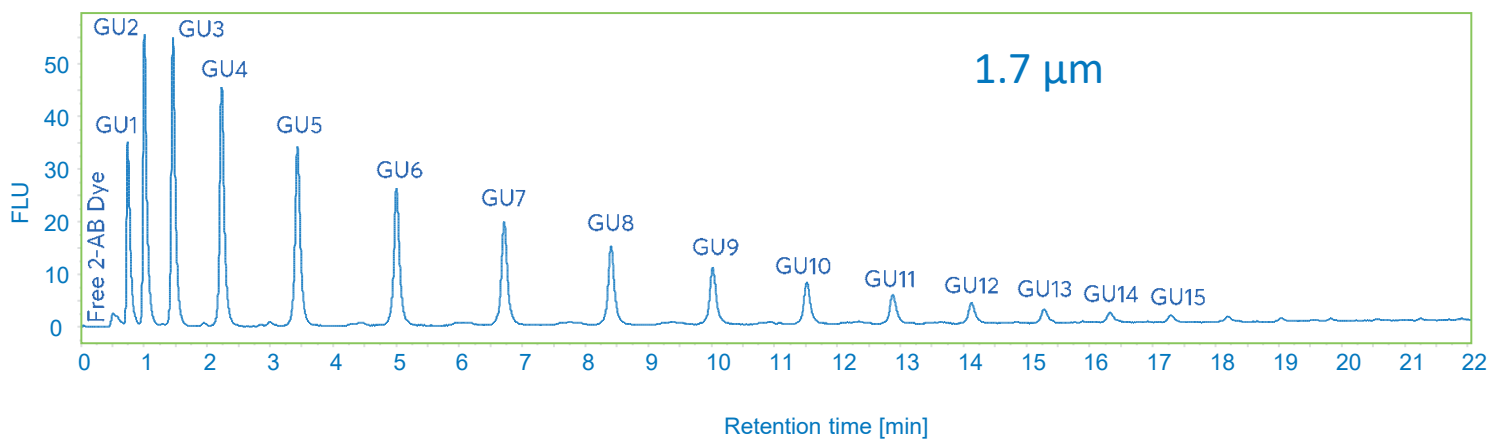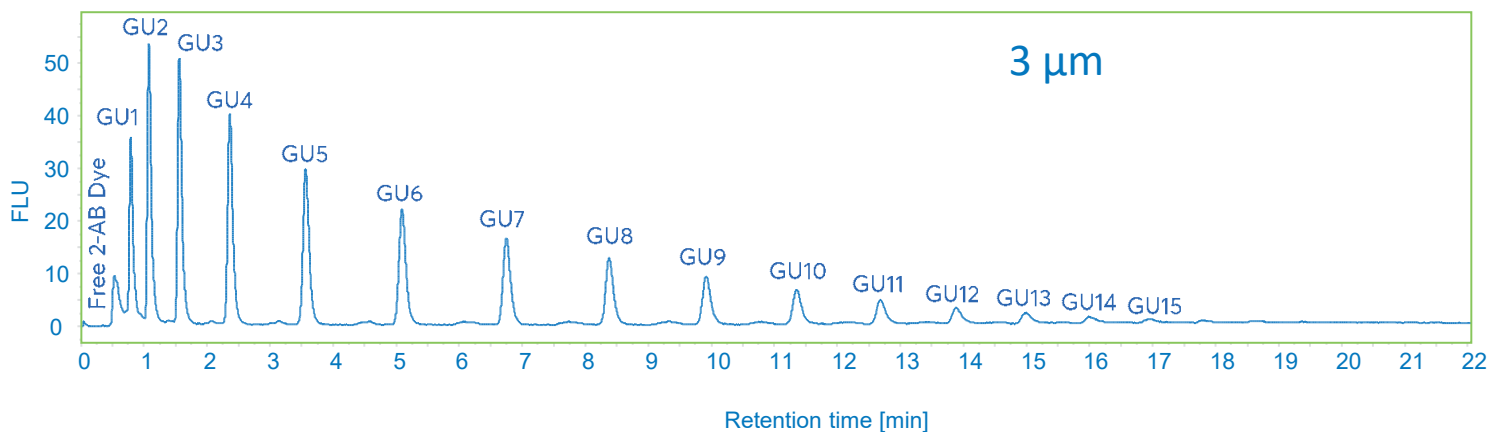


Sample: 2-AB labelled Dextran Ladder standard
Columns: ACE Excel 1.7 Glycan, 100x2.1 mm; ACE Excel 3 Glycan, 100x2.1 mm
Cat. No.: EXL-1716-1002 and EXL-1116-1002
Mobile phase: A = 100 mM ammonium formate (pH 4.5) and B = MeCN
Gradient: 75 – 52.5% B in 24 min
Flow rate: 0.5 ml/min
Temperature: 55 °C
Injection: 1 μl
Detection: FLD, λ_{ex} = 260 nm, λ_{em} = 430 nm

EXCELLENT COLUMN STABILITY

- The ACE® Excel® Glycan column shows good stability for glycan analysis over several thousand column volumes as seen in the graph below

COLUMN STABILITY UNDER GLYCAN ANALYSIS CONDITIONS

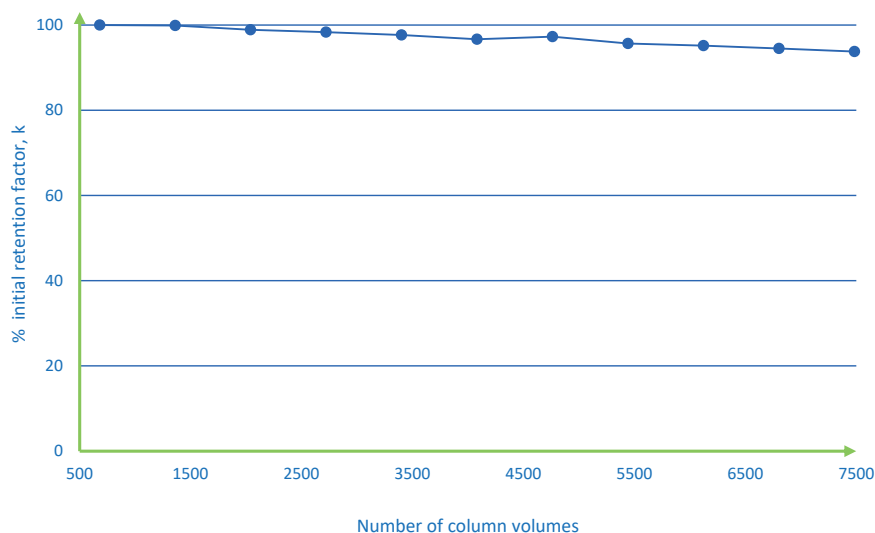

Sample: 2-AB labelled Dextran Ladder standard
Column: ACE Excel 1.7 Glycan, 50x2.1 mm
Cat. No.: EXL-1716-0502
Mobile phase: A = 100 mM ammonium formate (pH 4.5) and B = MeCN
Gradient: 75 – 60% B in 8 min
Flow rate: 0.5 ml/min
Temperature: 55 °C
Injection: 0.5 µl
Detection: FLD, λ_{ex} = 260 nm, λ_{em} = 430 nm

AVANTOR® ACE® EXCEL® GLYCAN UHPLC/HPLC ORDERING INFORMATION

Avantor® ACE® Excel® Glycan columns

Column dimensions	1.7 µm	3 µm
50 x 2.1 mm	EXL-1716-0502	EXL-1116-0502
100 x 2.1 mm	EXL-1716-1002	EXL-1116-1002
150 x 2.1 mm	-	EXL-1116-1502
50 x 4.6 mm	-	EXL-1116-0546
100 x 4.6 mm	-	EXL-1116-1046
150 x 4.6 mm	-	EXL-1116-1546

Avantor® ACE® Excel® Glycan guard cartridges

Description	Cat. No.
For analytical columns (2.1 mm ID)	EXL-1P16-GD2U
For analytical columns (4.6 mm ID)	EXL-1P16-GD4U

Related products

Description	Cat. No.
UHPLC guard holder	H0011
ACE Excel UHPLC column connector	EXL-CC10

Avantor® ACE®

CONTACT INFO:

vwr.com/ace
chromsupport@avantorsciences.com

About VWR, part of Avantor

As Avantor's channel brand, VWR serves as a leading global provider of product and service solutions to laboratory and production customers in the pharmaceutical, biotechnology, industrial, education, government and healthcare industries, and offers an integrated, seamless purchasing experience that is optimized for the way our customers do business. We set science in motion to create a better world. For more information, visit [avantorsciences.com](https://www.avantorsciences.com) or [vwr.com](https://www.vwr.com).



AUSTRIA

VWR International GmbH
Graumannsgasse 7
1150 Vienna
Tel.: +43 1 97 002 0
info.at@vwr.com

BELGIUM

VWR International bv
Researchpark Haasrode 2020
Geldenaaksebaan 464
3001 Leuven
Tel.: +32 (0) 16 385 011
vwr.be@vwr.com

CANADA

VWR International
2360 Argentia Road
Mississauga, Ontario L5N 5Z7
Tel.: +1 800 932 5000
Canada_Orders@vwr.com

CHINA

VWR (Shanghai) Co., Ltd
Bld.No.1, No.3728 Jinke Rd,
Pudong New District
Shanghai, 201203- China
Tel.: 400 821 8006
info_china@vwr.com

CZECH REPUBLIC

VWR International s. r. o.
Veetee Business Park
Pražská 442
CZ - 281 67 Stříbrná Skalice
Tel.: +420 321 570 321
info.cz@vwr.com

DENMARK

VWR International A/S
Tobaksvejen 21
2860 Søborg
Tel.: +45 43 86 87 88
info.dk@vwr.com

FINLAND

VWR International Oy
Valimotie 9
00380 Helsinki
Tel.: +358 (0) 9 80 45 51
info.fi@vwr.com

FRANCE

VWR International S.A.S.
Le Périgores – Bâtiment B
201, rue Carnot
94126 Fontenay-sous-Bois cedex
Tel.: 0 825 02 30 30* (national)
Tel.: +33 (0) 1 45 14 85 00 (international)
info.fr@vwr.com
* 0,18 € TTC/min + prix appel

GERMANY

VWR International GmbH
Hilpertstraße 20a
D - 64295 Darmstadt
Tel.: 0800 702 00 07* (national)
Tel.: +49 (0) 6151 3972 0 (international)
info.de@vwr.com
*Freecall

HUNGARY

VWR International Kft.
Simon László u. 4.
4034 Debrecen
Tel.: +36 52 521130
info.hu@vwr.com

INDIA
Avantor Performance Materials India Limited
17th Floor, Building No. 5, Tower C
DLF Cyber City Phase – III
Gurgaon - 122002, Haryana
Tel.: +91-1244-65-6700
chromatography.india@avantorsciences.com

IRELAND

VWR International Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 (0) 1 88 22 222
sales.ie@vwr.com

ITALY

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: +39 02 3320311
info.it@vwr.com

KOREA

VWR International ~
17, Daehak 4-ro, Yeongtong-gu,
Suwon-si, Gyeonggi-do
Tel.: +82 31 645 7256
saleskorea@avantorsciences.com

THE NETHERLANDS

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: +31 (0) 20 4808 400
info.nl@vwr.com

MEXICO

VWR International, S.de R.L. de C.V.
Km. 14.5 Carretera
Tlalnepantla-Cuautitlán
Col. Lechería
Tultitlán Edo. de México
CP 54940
Tel.: +52 (55) 5005 0100
vwrmx@vwr.com

MIDDLE EAST & AFRICA

VWR International FZ-LLC
Office 203, DSP Lab Complex,
Dubai Science Park,
Dubai, United Arab Emirates
Tel.: +971 4 5573271
Info.mea@vwr.com

NORWAY

VWR International AS
Brynsalleen 4,
0667 Oslo
Tel.: +47 22 90 00 00
info.no@vwr.com

POLAND

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: +48 58 32 38 200
info.pl@vwr.com

PORTUGAL

VWR International -
Material de Laboratório, Lda
Centro Empresarial de Alfragide
Rua da Indústria, nº 6
2610-088 Amadora
Tel.: +351 21 3600 770
info.pt@vwr.com

SINGAPORE

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel.: +65 6505 0760
sales.sg@vwr.com

SPAIN

VWR International Eurolab S.L.U.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: +34 902 222 897
info.es@vwr.com

SWEDEN

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: +46 (0) 8 621 34 00
kundservice.se@vwr.com

SWITZERLAND

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: +41 (0) 44 745 13 13
info.ch@vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard - Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: +44 (0) 800 22 33 44
uksales@vwr.com

UNITED STATES

VWR International, LLC
100 Matsonford Road
Building One Suite 200
Radnor, PA 19087
Tel.: +1 800 932 5000
VWRCustomerService@vwr.com

GO TO [VWR.COM](https://www.vwr.com) FOR THE LATEST NEWS, SPECIAL OFFERS AND DETAILS FROM YOUR LOCAL VWR SUPPORT TEAM

Prices, product, and/or services details are current when published and subject to change without notice. | Certain products or services may be limited by country, federal, state, provincial, or local regulations. | VWR, part of Avantor, makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC and/or Avantor, Inc. or affiliates. Offers valid in countries listed above, void where prohibited by law or company policy, while supplies last. | Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit [vwr.com](https://www.vwr.com) to view our privacy policy, trademark owners, and additional disclaimers.